

Product information

QF 24 V4
 V1 Jan 2018

Rapid Blood Genomic DNA Extraction Kit

Catalog #: Bt4782
Size: 50 preps
Storage: 4°C (1 year)

PRODUCT IS FOR SCIENTIFIC
 RESEARCH ONLY. NOT INTENDED
 FOR HUMAN OR ANIMAL USE.

Product Description:

The kit allows simple and fast isolation of high quality genomic DNA from fresh and anti-coagulated blood. The entire procedure takes approx. 15 minutes only. Up to 5 ml of blood sample can be treated for each mini-preps. Average DNA yields are 20~60 µg per ml of the whole blood sample. Purified DNA can be used in a wide range of downstream applications.

Contents:

Description	Size
Buffer TBP	40 ml
Universal Buffer Digestion	10 ml
Buffer PR	4 ml
Proteinase K	1.2 ml
TE Buffer	50 ml
Protocol	1

Features:

- **High quality:** OD260/OD280 of purified DNA is between 1.8~2.0, and can be directly used for downstream applications, such as PCR, Restriction Enzyme Digestion, Southern blotting, etc.
- **Simple procedure:** can be completed in 15 mins.
- **High Yield:** The yield of DNA is generally at 20~60 µg per 1 ml whole blood, and up to 5 ml of whole blood can be used in a single prep.
- Easy to scale up.
- Non toxic.

Protocol:

- Transfer 300µl anti-coagulated blood to a clean 1.5 ml tube. Add 600µl Buffer TBP, mix well and incubate for 1 min at room temperature.
Note 1: For frozen blood sample, pre-warm sample at 37°C for 3 minutes.
Note 2: For a fowl blood sample, transfer 100µl of the blood sample to 1.5 ml microtube, go to step 4.
- Centrifuge at 4000 x g for 1 min at room temperature, discard supernatant carefully.
- Add 500µl TE Buffer, suspend the deposit gently but thoroughly, Centrifuge at 4000 x g for 1 min at room temperature, discard supernatant carefully.
Note: Repeat this step once if necessary.
- Add 180µl Buffer Digestion and 20µl Proteinase K, mix thoroughly. Incubate at 56°C for 20~30 min.
Note: To obtain RNA-free DNA, add 20µl RNase A solution (20 mg/ml) to the tube, mix thoroughly and incubate at room temperature for 5 minutes.
- Add 60µl of Buffer PR, mix by inverting 5~10 times, incubate at -20°C for 5 minutes.
- Centrifuge at 12,000 x g for 5 minutes at room temperature. Transfer the supernatant to a new 1.5 ml tube.
- Add equal volume of isopropanol (approx 0.15~0.25 ml) to the solution, mix well by inverting 5 times. Incubate at room temperature for 2~5 minutes. Centrifuge at 12,000 x g for 5 minutes, discard the supernatant carefully.
- Add 1 ml of pre-cooled 75% ethanol to the tube, mix well by inverting 10 times. Centrifuge at 12,000 x g for 1 minute, discard the supernatant.
- Repeat the Step 8 once.
- Air-dry the pellet at room temperature with the lid open for 2~5 minutes.
Note: Don't over dry.
- Add 50~200µl of TE buffer to dissolve DNA pellet. Keep at 4°C for a couple hours until DNA pellet is completely dissolved. Purified DNA is ready for use. Or store at -20°C for long term storage.